

How to use the live microscope in the path conference room:

**Quick Start Guide to Live Microscope Display (DP70)**

1. Log into the Partners computer and turn on the Microscope. Make sure channel is set to “1” (on controller inside the podium) . Turn on the projector with the remote (RGB1).
2. On the computer desktop find and double click the ‘live Micro icon’. The microscope viewer software will open. (Olympus DP Controller).
3. **( F7) In the viewing software Press the F7 key on the keyboard this will start the live mode.**
4. **White balance. The white balance eye dropper button is on the top row of menu icons. Click white balance button and then with the tool click on a white or clear area.**
5. **(F11) Press F11 key on the key board the active the Live Full Screen mode. This is the mode used to view your sides.**
6. When done with the viewing software press the F11 or Esc button and exit the software.
7. Turn off the microscope and computer when done.